

QUANTOM™ Utility

Product Description

QUANTOM™ Utility software is used to access data saved to the QUANTOM Tx™ Microbial Cell Counter from a remote PC via WiFi.

PC Requirements

OS: Windows 10, 32-bit
Display: 1280 x 1024 pixels
Internet: WiFi connection

Directions for Use

CONNECT THE QUANTOM Tx™ TO WIFI


1. Connect the supplied WiFi dongle to a USB port.
2. Open WiFi Options. Log in to a wireless network.
3. Make note of the IP address.

INSTALL QUANTOM™ UTILITY ON A PC

1. Download QUANTOM™ Utility from www.logosbio.com.
2. Unzip the file and install on a compatible PC.

CONNECT QUANTOM™ UTILITY TO THE QUANTOM Tx™


1. Open QUANTOM™ Utility.

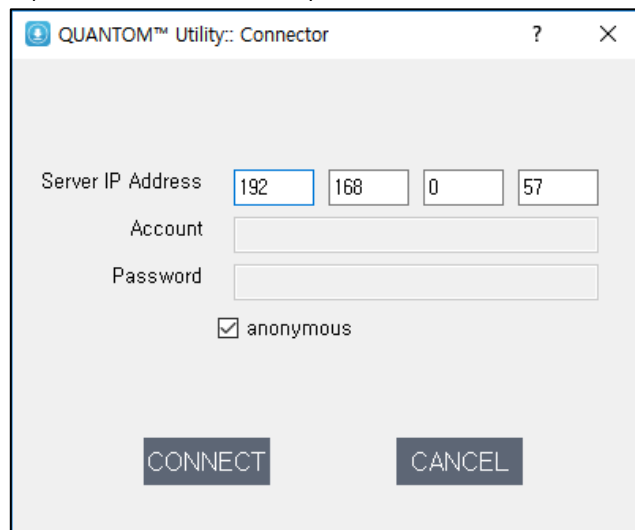

2. Enter the QUANTOM Tx™ IP address. Press **CONNECT**.

- ! The QUANTOM Tx™ and PC must be connected to the same wireless network.
- ! If the 21 CFR Part 11 Package (Q11001) has been installed to the QUANTOM Tx™, a user name and password will need to be entered to access data.

3. Select folders/files saved to the QUANTOM Tx™.

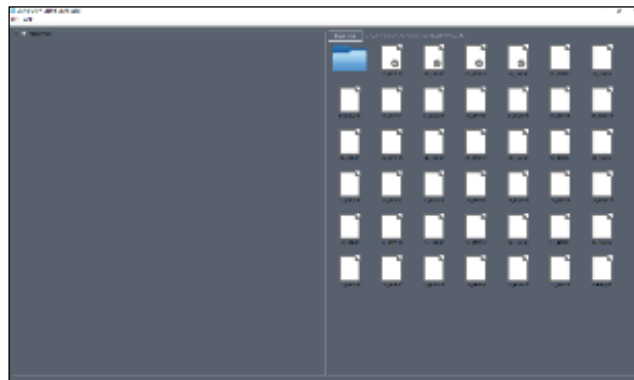

4. Press **Download** to import to your PC.



HEADQUARTERS

Doosan Venturedigm Ste 514
415 Heungan-daero, Dongan-gu
Anyang-si, Gyeonggi-do 14059
South Korea

Tel: +82 (31) 478-4185

USA

7700 Little River Turnpike Ste 207
Annandale, VA 22003
USA

Tel: +1 (703) 622-4660

EUROPE

11B avenue de l'Harmonie
59650 Villeneuve d'Ascq
France

Tel: +33 (0)3 74 09 44 35