

Electrophoretic Tissue Clearing Solution

C13001

Storage

Room temperature

Product Description

Appearance	Colorless liquid
pH	8.5

Electrophoretic Tissue Clearing Solution is a premixed SDS-based buffer optimized for use with the X-CLARITY™ Tissue Clearing System.

Directions for Use

If precipitates form, incubate at 60°C for 1 hour or at room temperature for 1 day.

1. Use 1.2 L fresh Electrophoretic Tissue Clearing Solution for each run with the X-CLARITY™ Tissue Clearing System.
2. Replace the solution every 24 hours.

! *Reused clearing solution may have lowered buffering capacity, which reduces clearing efficiency. Prolonged use of Electrophoretic Tissue Clearing Solution may cause electrolysis of the extracted lipids, turning the tissue yellow.*

Disclaimer

This product is for research use only. Please consult the material safety data sheet for information regarding hazards and safe handling practices.

References

1. Lee, E et al. ACT-PRESTO: Rapid and consistent tissue clearing and labeling method for 3 dimensional (3D) imaging. *Sci Rep* 6, 18631 (2016).
2. Yang, B. et al. Single-cell phenotyping within transparent intact tissue through whole-body clearing. *Cell* 158, 945–958 (2014).
3. Chung, K. et al. Structural and molecular interrogation of intact biological systems. *Nature* 497, 332–337 (2013).

Additional information is available on our website at www.logosbio.com.
© Logos Biosystems 2018. All rights reserved.



biosystems

HEADQUARTERS

FL 2 & 3
28 Simindaero 327beon-gil, Dongan-gu
Anyang-si, Gyeonggi-do 14055
South Korea

Tel: +82 (31) 478-4185

USA

7700 Little River Turnpike STE 207
Annandale, VA 22003
USA

Tel: +1 (703) 622-4660, +1 (703) 942-8867

EUROPE

11B avenue de l'Harmonie
59650 Villeneuve d'Ascq
France

Tel: +33 (0)3 74 09 44 35