


This Installation Guide provides instructions for installing the CELENA<sup>®</sup> S Digital Imaging System. For detailed instructions, refer to the user manual in the supplied USB drive.

### PLACEMENT


Lift the CELENA<sup>®</sup> S out of its box by grasping its base firmly with both hands at diagonally opposite ends. Place on a clean, level, and sturdy surface.

- ! Do not hold or lift the CELENA<sup>®</sup> S by its top cover.
- ! Avoid vibrations from other devices.
- ! Leave sufficient space around the instrument for proper ventilation and to prevent overheating.
- ! Do not expose the instrument to intense ultraviolet light.


### STAGE ASSEMBLY

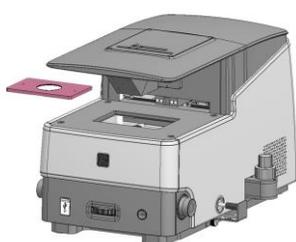

1. Place the insert plate onto the stage with the flat side up.

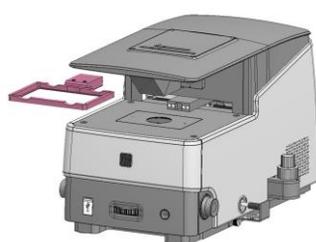


2. Attach the vessel holder frame to the vessel holder frame arm. The frame will snap into place.

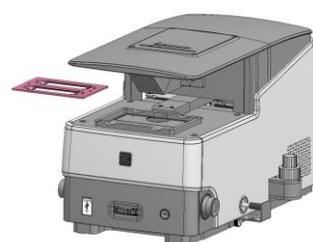


3. Place the desired vessel holder into the vessel holder frame.

### SELECTION TAIL GUARD

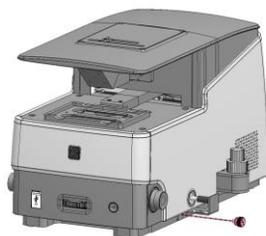


1. Unscrew the screw from the selection tail guard.

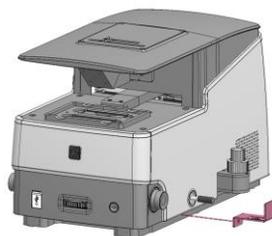

2. Pull the guard away from the instrument. Store in a safe place for future use.

### CONNECTIONS

Make the following connections at the back of the CELENA<sup>®</sup> S.

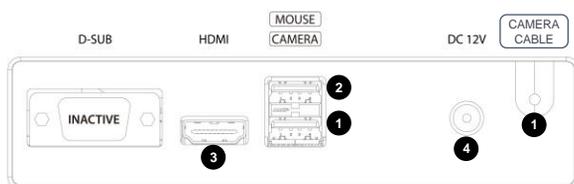

1. Insert the camera cable into the *bottom* USB port.
2. Plug the wireless mouse dongle into the *top* USB port.
3. Connect an HDMI-compatible monitor to the CELENA<sup>®</sup> S via the HDMI cable.
4. Connect the power cord to the DC 12V power inlet and an electrical outlet.

\*For the EU version of the CELENA<sup>®</sup> S, the camera cable is internalized and the inactive D-SUB port has been removed.