


CountWire™

Version Release Note: 1.3.6

CountWire

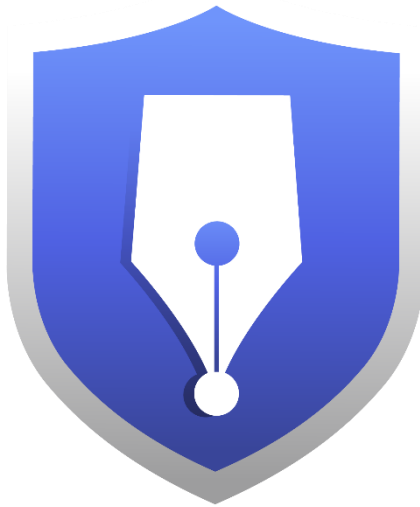

Jun. 11, 2025

Sales & Marketing Department

< Version 1.3.6 >

Release Date	Model	Version	Writer
Jun. 11, 2025	CountWire™ Client	1.3.6	YuBeen Lee

Hardware	Software	New features	Improvement
-	o	-	o

Version description

<Bug fix>

1. Disabled FL-related settings when BX7 model is selected

- Adjusted the configuration so that all FL-related options are automatically disabled when the BX7 model is selected.

2. Reset user password

- Addressed an issue where resetting a user's password would fail if the minimum length requirement was set between 17 and 20 characters.
- Improved the password reset feature to create a random password that follows the minimum length setting, instead of always using a fixed 8-character password.

<Update>

3. Spectra Max slide

- Added a Spectra Max option to the available slide types.
- The Spectra Max Option utilizes the same slide as the Spectra Option. The distinction between the two lies in the extent of the imaging area.

4. Strengthened password generation logic

- Modified the auto-generation algorithm to enforce inclusion of all required character types—uppercase, lowercase, numeric, and special—for improved security.

5. Improved Slide option behavior based on connected LUNA-BX7/FX7 model SW version

- If the connected software version is higher than 1.9.5: Spectra Max and Spectra options are enabled.
- If the version is exactly 1.9.5: Spectra Max is disabled and Spectra is enabled.
- If the version is lower than 1.9.5: both Spectra Max and Spectra are disabled.

6. Log event when Delete action is performed

- Added functionality to record an event log entry whenever the Delete action is executed from the Report list.

Please note!

To ensure proper compatibility, please use the following versions:

- LUNA-FX7: version 1.9.6
- CountWire Client: version 1.3.6
- CountWire DSS: version 1.3.5

*Download the latest software via website <https://logosbio.com/support/software>

- End of Page -