

CountWire™

Version Release Note: 1.3.5

CountWire



Apr. 02, 2025

Sales & Marketing Department

< Version 1.3.5 >

Release Date	Model	Version	Writer
Apr. 02, 2025	CountWire™ Data Storage Server	1.3.5	YuBeen Lee

Hardware	Software	New features	Improvement
-	o	-	o

Version description

<Bug fix>

1. Protocol

- Addressed an issue where excessively long protocol names were not fully displayed in the list view.
- Standardized the maximum protocol name length to 37 characters, aligning with the specification used in LUNA-FX7.
- Enhanced the protocol overwrite validation to include comparisons with existing items in the protocol list.
- Implemented a feature to check whether a protocol has been modified by another client prior to saving.

<Update>

2. Spectra slide

- Added a Spectra option to the available slide types.
- Introduced a dedicated protocol for spectra, SPECTRA_AP-1, available in FL mode.

3. Online installer(.dss)

- The installer has been converted to an online installer, resulting in a reduced file size compared to the previous version.
- Mitigated potential compatibility issues that may arise when running the software on newer versions of Ubuntu.

4. Time sync

- Added a notification popup to alert users when there is a time mismatch between the client and the server systems.

5. User Interface Enhancements

- Optimized the performance of the Report List display for faster loading times.
- Improved the user interface to improve overall usability.

Please note!

To ensure proper compatibility, please use the following versions:

- LUNA-FX7: version 1.9.6
- CountWire Client: version 1.3.6
- CountWire DSS: version 1.3.5

*Download the latest software via website <https://logosbio.com/support/software>

- End of Page -