

# **LUNA-FX7™ Automated Cell Counter**

**Version Release Note: 1.9.6**



**Jun. 19, 2025**

**Sales & Marketing Department**

## < Version 1.9.6 >

Release Date	Model	Version	Writer
Jun. 28, 2025	LUNA-FX7™	1.9.6	Daegeun Han
Hardware	Software	New features	Improvement
-	o	-	o

### Version description

#### <Bug fix>

##### 1. Print

- Fixed an issue where buttons could be pressed when the printer device was not connected.
- Fixed an issue where Reusable/Validated/Spectra slide types would appear as 9/10/11 channel slides.
- This bug has been fixed.

#### <Update>

##### 1. Spectra slide

- Change the insertion position of the slide: Move the existing FOV forward by one space.
- Capture and count 5 images in total, FOV range 6~10 of Spectra slide.

##### 2. Spectra Max slide

- Spectra Max slide option added.
- Like previous versions of Spectra slide, it acquires and counts 10 images in total from 1 to 10.
- 

##### 3. Changes the exponent notation for Counting results.

- $X < 1.9.6$  :  $8.93 \times 10^6$  cells/mL
- $1.9.5 < X$  :  $8.93 \times 10^6$  cells/mL

#### **Note for CountWire customers!**

To ensure proper compatibility, please update CountWire to version 1.3.6 when updating LUNA-FX7 to version 1.9.6.

Download the latest software via website <https://logosbio.com/support/software>

**- End of Page -**