

# **LUNA-FX7™ Automated Cell Counter**

**Version Release Note: 1.9.5**



**Mar. 28, 2025**

**Sales & Marketing Department**

## < Version 1.9.5 >

Release Date	Model	Version	Writer
Mar. 28, 2025	LUNA-FX7™	1.9.5	Yubeen Lee

Hardware	Software	New features	Improvement
-	o	-	o

### Version description

#### <Bug fix>

1. Counting Result(CSV file)
  - An issue occurred where the PDF changed after approval due to a missing title in the CSV file.
  - This bug has been fixed.
2. Auto logout
  - An issue occurred where the system would automatically log out during tasks such as counting and saving.
  - This bug has been fixed.

#### <Update>

1. Spectra slide
  - A spectra option has been added to the slide options.
  - A protocol for spectra (SPECTRA\_AP-1) has been added in FL mode..
2. Safe position
  - To prevent the slide from falling, a new feature has been added that moves the slide insertion part to a safe position if it remains idle for a certain period...

Download the latest software via website <https://logosbio.com/support/software>

**- End of Page -**