

# QUANTOM Tx™ Microbial Cell Counter

**Version Release Note: 1.9.2**



**Oct. 7, 2021**

**Sales & Marketing Department**

## < Version 1.9.2 >


Release Date	Model	Version	Writer
Oct. 7, 2021	QUANTOM Tx™	1.9.2	Deokhui Chang

Hardware	Software	New features	Improvement
-	0	-	0

### Version description

1. New WiFi dongle is supported
  - Network connection of old type WiFi dongle was reported to be unstable.
  - WiFi dongle was changed to new type and related software has been updated.
2. Roundness calculation formula was modified
  - Roundness calculation formula was changed for more accurate calculation of rod-shaped bacterial cells
3. Improved calibration routine
  - Fixed the problem of failing when performing calibration of some under-powered LEDs.
  - When calibration is complete, the calibrated exposure value is displayed in the calibration values.

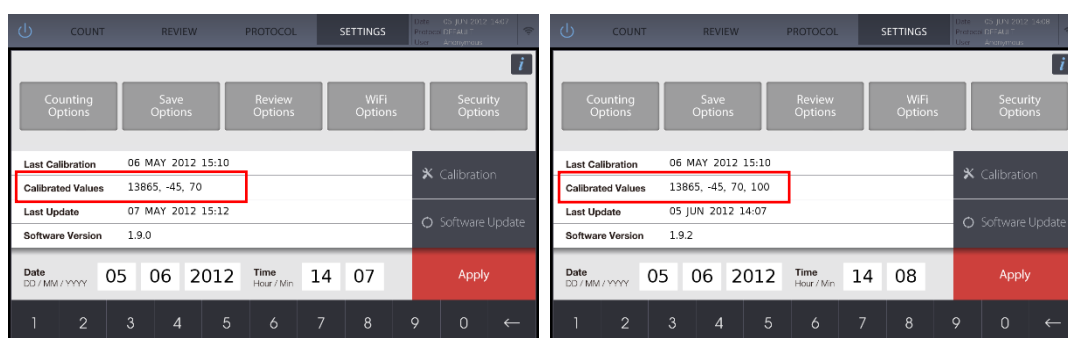

**⚠ Do not update if the current version is 1.9.3 or 1.9.4.**

- End of Page -